

Tissue Processor MTP



... the most universal unit
easy to operate

The tissue processor program meets all the requirements of today's histo-pathology, covering safety of the specimen, fast processing, flexibility, environmental protection and economy,

9001: 2000

MTP – specification

Protection hood: serves for covering moveable equipment parts during operation and as suction hood for stopping bad odours, a side opening in the hood for container loading

Glass containers: 2 litres with holder as well as a stop peg.

Steel container: 2 litre contents, heated. double walled, heating capacity amounts to 70 W with 24 volts DC, temperature measured by sensor and for overtemperature protection a safety temperature limitation is built in, temperature value can be set for every heated container separately.

Stainless steel baskets: for receiving the tissue samples, filled outside, closed with a snap lock, filled baskets can be loaded to the receiving arms

Lifting unit: lifting and lowering loaded baskets and closing lids available for each container

Rotation unit: moves all containers in the programmed sequence

Control unit: LCD display with 4 lines and only 5 operating keys, free programmable, possibility to store up to 10 programs with individual rest and dripping time for all baths, present active baths, remaining rest time, possibility of a start delay of up to 99 hours

Extraction : rear connection possibility for a suction hose, 100 mm diam, optionally the unit can be equipped with its own exhaust fan as well as charcoal filter.

Electrical data: 100 – 265 V / 50/60Hz

Dimensions/
Weight: 87 x 87 x 70 cm - approx. 65 kg



SLEE

MAINZ

SLEE MEDICAL GMBH
P. O. Box 10 05 07
D-55136 Mainz – Germany
FON (+49) 6131/9 58 71-0
FAX (+49) 6131/9 58 71-22

Your authorised distributor

ML chemica
Zahradní 1a
664 41 Troubsko u Brna

tel.: 547 227 533
tel., fax: 547 227 532

e-mail: info@mlchemica.cz

www.mikroskopy.net